FRANKA HAND

QUICK GUIDE FOR INSTALLATION

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NOTICE

This guide is not the manual of your Franka Hand and does not replace it.

INDICATIONS

Before operating the Hand, all persons working with it have to read the corresponding manual carefully. Read or download the manual, also in other languages, here: www.franka.de/documents.

Don't hesitate to contact your sales partner or integrator if you need more detailed information. Alternatively, you can get in touch with us directly by sending an e-mail to support@franka.de.

NOTICE

Please carefully follow all instructions, warnings and other notes in this document and the official manual. Any deviation from the steps prescribed in these documents constitutes a misuse (as described in the paragraph "Misuse" of the corresponding manual). Franka Robotics does not assume any warranty and/or liability for any possible consequences.

Get your manual and additional supporting material
in English and other languages.
www.franka.de/documents

CONNECTING THE HAND

**This booklet will guide you
through the product manual's steps
to connect the Hand.**

Document number: R50040
Release Version: 1.1

The English documentation is the ORIGINAL DOCUMENTATION.
Other languages are translations of the original document.

See Franka Hand product manual
Paragraph: About Franka Hand

NOTICE

Always keep the original packaging in case of relocating the Hand.

NOTICE

Franka Hand is not part of the certified machinery.

About Franka Hand

1	Finger
2	Fingertip
3	Plug
4	Screw holes
5	Flange to Arm
6	Hole for alignment pin

Franka Hand is an electrical two-finger parallel gripper produced by Franka Robotics. No external cabling is needed. Communication and power supply for the Hand is established via the flange connector of the Arm.

Main components:

- Franka Hand

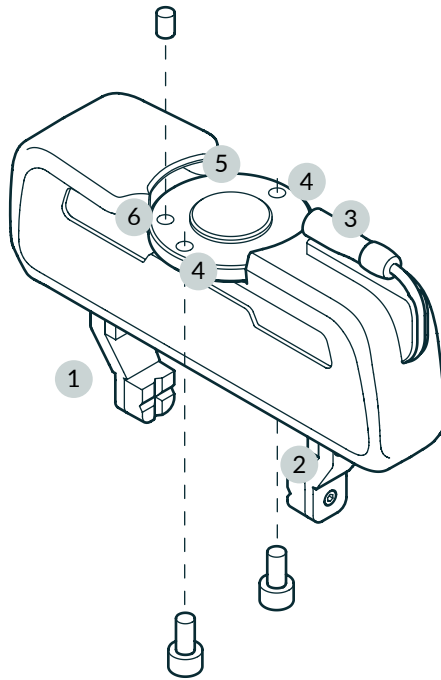
Also included:

- 1 set of fingertips
- 2x DIN7984 M6x12 ST 8.8 screws
- 1x ISO2338B 6x10 h8 A2 cylindrical pin

Required equipment (*not included*):

- Hex key size 4
- A robot system of Franka Robotics

Overview

See Franka Hand product manual
Paragraph: Mounting the Hand

⚠ WARNING!

Falling Hand

Risks of severe injury, such as crushing to fingers, hands, upper body parts, head.

- Ensure the Hand is mounted correctly.

⚠ WARNING!

Risk of electric shock

Risk of personal injury by electric shock.

- Disconnect the Arm from the mains power supply before mounting the Hand.

⚠ WARNING!

Falling and/or flying tools from the robot

Tools that remain inserted in the robot may turn into projectiles during later motions of the Arm and lead to injuries.

- Do not leave any tools inside the robot.

Mounting and connecting the Hand

Required material:

- 2x M6x12 screws
- 1x cylindrical pin

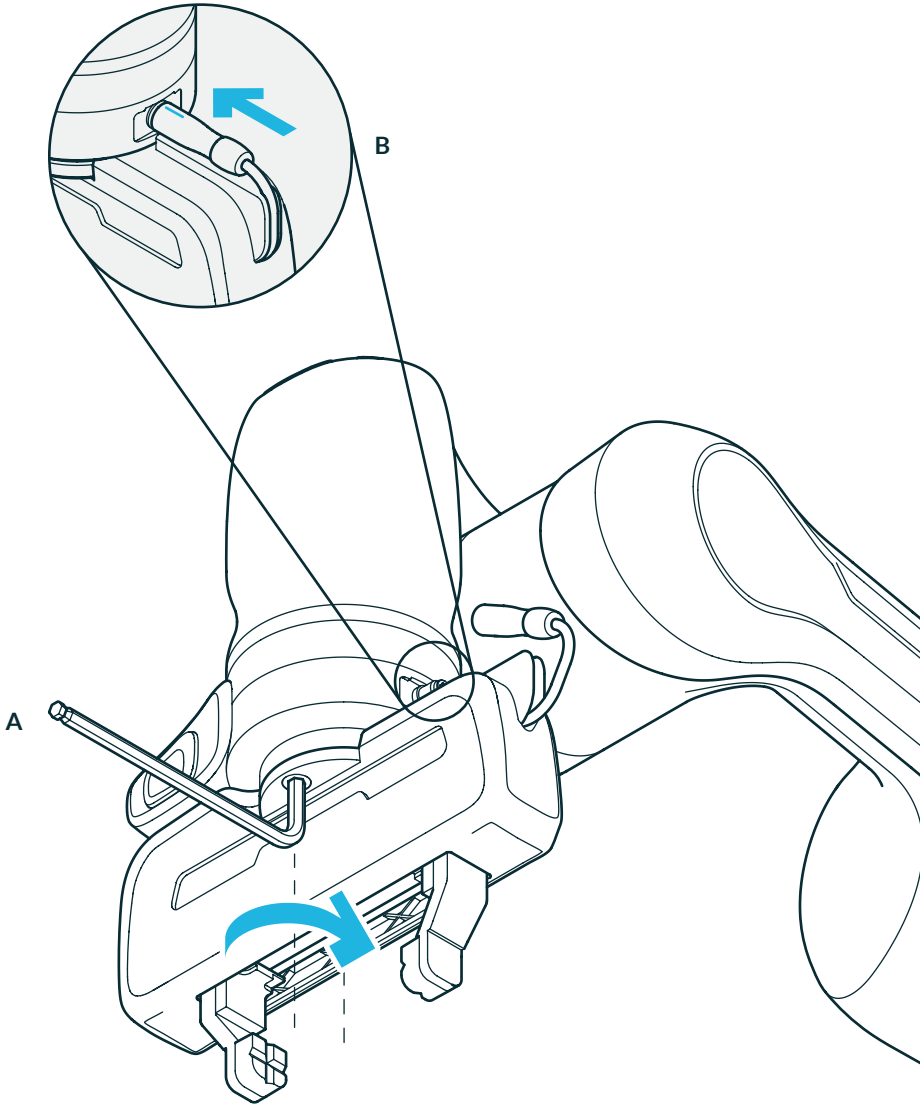
Precondition:

- The robot is switched off

Procedure:

1. If desired, insert the cylindrical pin in the H7 fit of the flange of the Hand.
2. Position the Hand with its flange on the end effector flange on the Arm. Should you not use the cylindrical pin, make sure that the side of the Hand with the cable is attached on the side of the Pilot-Grip where the connector is located.
3. Attach the Hand with 2x M6x12 screws (do not use other types of screws) and use 5 Nm of fastening torque (A).
4. Insert the plug connector of the Hand into the X6 connector on the Pilot-Grip of the Arm (B).

Step 1

See Franka Hand product manual
Paragraph: Configuration

WARNING!

Unexpected motions during guiding due to incorrectly configured end effectors

Incorrectly configured end effector mass and inertia may result in the gravitational forces not being fully compensated for. The resulting and unexpected behaviour of the robot including end effector may lead to injuries such as crushing, tearing of skin, and puncturing.

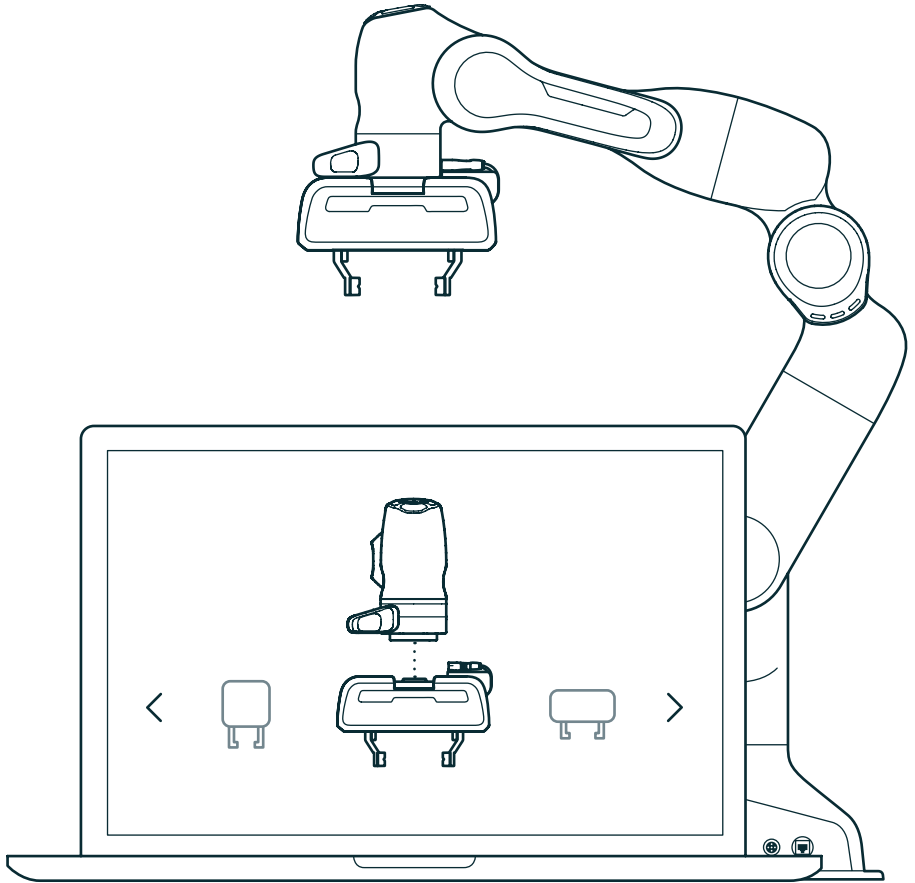
- Always check the configuration of the end effector.
- When copying an already parametrized app or task to another Franka Robotics Production system, ensure that the end effector configuration is still identical to the original one.

Configuring the Hand

When first starting a Franka robot - by means of the initial configuration - or at a later point in time by accessing the end effector settings, the Hand can be selected from a dedicated drop down menu.

For further information please refer to the relevant robot's product manual of Franka Robotics.

Step 2

See Franka Hand product manual
Paragraph: Using the Hand

WARNING!

Falling objects from end effectors due to loose coupling or malfunction

Objects falling from the gripper leads to injuries to hands, fingers, feet, and toes.

- Always wear personal protective equipment (e.g., safety shoes).
- Use the appropriate type of grippers to prevent objects from falling.
- Consider the shape, texture, and weight of the grasped objects in the risk assessment according to ISO 10218-2.
- Using light and/or round objects may reduce risks significantly.

INDICATIONS

If you need more detailed information, please contact your sales partner or integrator. Alternatively you can get in touch with us directly by sending an e-mail to support@franka.de.

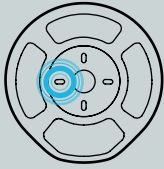
Using the Hand

By setting the Pilot-Mode button to control the end effector, the fingers of the Hand can be steered via the arrow keys of the Pilot-Disc:



- To slowly open the Hand, keep the “left” key pressed (A).
- To slowly close the Hand, keep the “right” key pressed (B).
- To open the Hand to full width, press the “down” key once (C).
- To make the Hand grasp with its holding force, press the “up” key once (D).

Quick tips


A



B

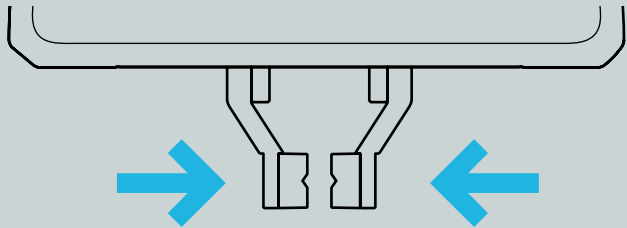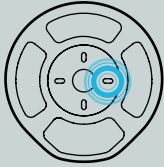C

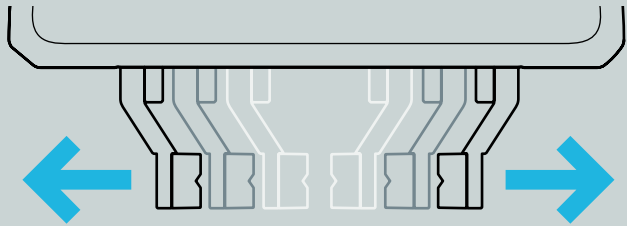D

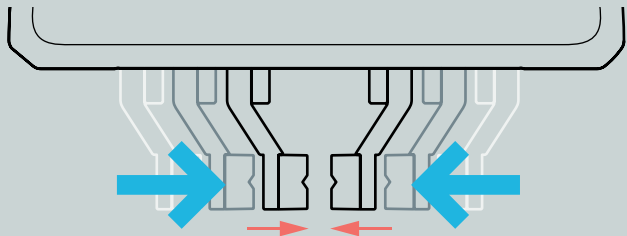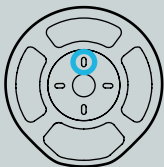Enjoy using your Franka Robotics robot. We look forward to getting to know about your success! Share your experience with us as a post, and we will amplify it on our own digital and social media platforms.

Just tag **@frankarobotics** or **#frankarobotics** in your post, and we will make sure your endeavors will get international visibility!

For more marketing and PR opportunities, reach out to **marketing@franka.de**.

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Franka Robotics GmbH